
pylinkwrapper Documentation

Release 1.0

Nick DiQuattro

March 16, 2016

1	Installation	3
2	Connector Class	5
3	Sample Experiment	9
4	Indices and tables	11
	Python Module Index	13

Contents:

Installation

1. [Download](#) and place the pylinkwrapper folder somewhere you can easily reach in the file system.
2. Add the folder to Psychopy's python path as detailed here: <http://www.psychopy.org/recipes/addCustomModules.html>
3. Update pylink to the latest version provided by the SR Research development kit. It should be installed on the display computer.

That's it! Here's a quick demo of getting connected to the Eyelink:

```
import pylinkwrapper

win = visual.window(monitor='nickMon', fullScr=True, allowGUI=False, color=-1)
tracker = pylinkwrapper.connect(win, '1_nd')
```

Check out the documentation for the [Connector Class](#) class to see what functions are available.

Connector Class

This class provides methods for interacting with the EyeLink from psychopy.

class `connector.Connect` (*window*, *edfname*)

Provides functions for interacting with the EyeLink via Pylink.

Parameters

- **window** – Psychopy window object.
- **edfname** (*str*) – Desired name of the EDF file.

calibrate (*cnum=13*, *paval=1000*)

Calibrates eye-tracker using psychopy stimuli.

Parameters

- **cnum** (*int*) – Number of points to use for calibration. Options are 3, 5, 9, 13.
- **paval** (*int*) – Pacing of calibration, i.e. how long you have to fixate each target in milliseconds.

convert_coords (*x*, *y*, *to='eyelink'*)

Converts from degrees visual angle units to EyeLink Pixel units.

Parameters

- **x** (*float or int*) – X coordinate in visual angle.
- **y** (*float or int*) – Y coordinate in visual angle.
- **to** – Direction of conversion. Options: 'eyelink' or 'psychopy'.

Returns Two values in order x, y

draw_ia (*x*, *y*, *size*, *index*, *color*, *name*)

Draws square interest area in EDF and a corresponding filled box on eye-tracker display. Must be called after `set_trialid()` for interest areas to appear in the EDF.

Parameters

- **x** (*float or int*) – X coordinate in degrees visual angle for center of check area.
- **y** (*float or int*) – Y coordinate in degrees visual angle for center of check area.
- **size** (*float or int*) – length of one edge of square in degrees visual angle.
- **index** (*int*) – number to assign interest area in EDF
- **color** (*int*) – color of box drawn on eye-tracker display (0 - 15)

- **name** (*str*) – Name of interest area in EDF

draw_text (*msg*)

Draws text on eye-tracker screen.

Parameters **msg** (*str*) – Text to draw.

end_experiment (*spath*)

Closes and transfers the EDF file.

Parameters **spath** (*str*) – Absolute file path of where to save the EDF file.

fix_check (*size, ftime, button*)

Checks that fixation is maintained for certain time.

Parameters

- **size** (*float or int*) – Length of one side of box in degrees visual angle.
- **ftime** (*float*) – Length of time to check for fixation in seconds.
- **button** (*char*) – Key to press to recalibrate eye-tracker.

get_gaze ()

Gets current gaze position of eye. Must be called between `record_on()` and `record_off()`. Sendlink must be set to True as well.

Returns list of coordinates in the form of [x, y].

record_off ()

Stops recording.

record_on (*sendlink=False*)

Starts recording. Waits 50ms to allow eyelink to prepare.

Parameters **sendlink** (*bool*) – Toggle for sending eye data over the link to the display computer during recording.

sac_detect (*x, y, radius*)

Checks if current gaze position is outside a circular interest area.

Parameters

- **x** (*float or int*) – X coordinate in degrees visual angle for center of circle IA.
- **y** (*float or int*) – Y coordinate in degrees visual angle for center of circle IA.
- **radius** (*float or int*) – Radius of detection circle in degrees visual angle.

Returns List of whether saccade was detected and the gaze coordinates

at time of detection. In the form of [bool, [x, y]]. :rtype: bool, list

send_command (*cmd*)

Sends a command to the Eyelink.

Parameters **cmd** (*str*) – Command to send.

send_message (*txt*)

Sends a message to the tracker that is recorded in the EDF.

Parameters **txt** (*str*) – Message to send.

send_var (*name, value*)

Sends a trial variable to the EDF file.

Parameters

- **name** (*str*) – Name of variable.
- **value** (*float, str, or int*) – Value of variable.

set_status (*message*)

Sets status message to appear while recording.

Parameters **message** (*str*) – Text object to send, must be < 80 char

set_trialid (*idval=1*)

Sends message that indicates start of trial in EDF.

Parameters **idval** – Values to set TRIALID.

set_trialresult (*rval=0, scrcol=0*)

Sends trial result to indicate trial end in EDF and clears screen on EyeLink Display.

Parameters

- **rval** (*float, str, or int*) – Value to set for TRIAL_RESULT.
- **scrcol** (*int*) – Color to clear screen to. Defaults to black.

Sample Experiment

This is a very simple experiment that demonstrates use of `pylinkwrapper`. It can be found in the `sample` folder.

```
"""
Pylink Wrapper test experiment
N. DiQuattro - January 2015

This is a simple experiment where a circle appears randomly on the screen. It's
purpose is to provide examples of how to use the pylink wrapper with a psychopy
experiment.

There's help documentaiton available for each of the functions that show the
available parameters.
"""

# Import modules
from psychopy import visual
from psychopy import core, event
import numpy as np

import pylinkwrapper # Here's the special one

# Window set-up
win = visual.Window(monitor='nickMon', units='deg', fullscr=True,
                    allowGUI=False, color=0)

# Initiate eye-tracker link and open EDF
tracker = pylinkwrapper.Connect(win, '1_test')

# Calibrate eye-tracker
tracker.calibrate()

# Stimulus
fix = visual.Circle(win, radius=1, pos=(0, 0), fillColor=[1, 0, 0],
                    lineColor=[1, 0, 0])

cfix = visual.Circle(win, radius=.15, fillColor=-1, lineColor=-1)

# Display stimulus 5 times
for t in range(5):
    # Find random coordinates and set them
    fx = np.random.randint(-10, 10)
    fy = np.random.randint(-10, 10)
    fix.setPos((fx, fy))
```

```
# Eye tracker trial set-up
stxt = 'Trial %d' % t
tracker.set_status(stxt) # Define status message that appears on eye-link
# display

tracker.set_trialid() # Sends trial start message for EDF
tracker.send_message('Circle Trial')

# Draw IA
tracker.draw_ia(fx, fy, 1, 1, 5, 'circle') # Draw interest area and box

# Start recording
tracker.record_on()

# Draw and display circle
cfix.draw()
fix.draw()
win.flip()

# Wait for response
keyp = event.waitKeys()

# Stop Recording
tracker.record_off()

# Send response key to EDF file
tracker.send_var('response', keyp[0][0])

# End trial for EDF
tracker.set_trialresult()

# ISI with fixation check
cfix.draw()
win.flip()
tracker.fixcheck(2, 1, 'z')

# Retrieve EDF
tracker.end_experiment('C:\\edfs\\') # Closes and retrieves EDF file to
# specified path
```

Indices and tables

- `genindex`
- `modindex`
- `search`

C

connector, 5

C

calibrate() (connector.Connect method), 5
Connect (class in connector), 5
connector (module), 5
convert_coords() (connector.Connect method), 5

D

draw_ia() (connector.Connect method), 5
draw_text() (connector.Connect method), 6

E

end_experiment() (connector.Connect method), 6

F

fix_check() (connector.Connect method), 6

G

get_gaze() (connector.Connect method), 6

R

record_off() (connector.Connect method), 6
record_on() (connector.Connect method), 6

S

sac_detect() (connector.Connect method), 6
send_command() (connector.Connect method), 6
send_message() (connector.Connect method), 6
send_var() (connector.Connect method), 6
set_status() (connector.Connect method), 7
set_trialid() (connector.Connect method), 7
set_trialresult() (connector.Connect method), 7