
phconvert Documentation

Release 0.7.2+0.g42d7ece

Antonino Ingargiola

Apr 13, 2017

Contents

1	Module <code>hdf5</code>	3
2	Module <code>loader</code>	7
3	Module <code>pqreader</code>	9
4	Module <code>bhreader</code>	11
5	Indices and tables	15
	Python Module Index	17

Version 0.7.2+0.g42d7ece ([release notes](#))

`phconvert` is a python 2 & 3 library which helps writing valid `Photon-HDF5` files. This document contains the API documentation for `phconvert`.

The `phconvert` library contains two main modules: *hdf5* and *loader*. The former contains functions to save and validate `Photon-HDF5` files. The latter, contains functions to load other formats to be converted to `Photon-HDF5`.

The `phconvert` repository contains a set the notebooks to convert existing formats to `Photon-HDF5` or to write `Photon-HDF5` from scratch:

- [phconvert notebooks \(read online\)](#).

In particular see notebook [Writing Photon-HDF5 files \(read online\)](#) as an example of writing `Photon-HDF5` files from scratch.

Finally, `phconvert` repository contains a [JSON specification](#) of the `Photon-HDF5` format which lists all the valid field names and corresponding data types and descriptions.

Contents:

The module `hdf5` defines functions to save and validate Photon-HDF5 files. The main two functions in this module are:

- `save_photon_hdf5()` to save data from a dictionary to Photon-HDF5.
- `assert_valid_photon_hdf5()` to validate if a HDF5 file is valid Photon-HDF5.

This module also provides functions to save free-form dict to HDF5 (`dict_to_group()`) and read a HDF5 group into a dict (`dict_from_group()`). Finally there are utility functions to easily print HDF5 nodes and attributes (`print_children()`, `print_attrs()`).

For more info see: [Writing Photon-HDF5 files](#).

List of functions

Main functions to save and validate Photon-HDF5 files.

```
phconvert.hdf5.save_photon_hdf5(data_dict, h5_fname=None, user_descr=None,
                                write=False, compression={'complib': 'zlib', 'com-
                                plevel': 6}, close=True, validate=True, warnings=True,
                                skip_measurement_specs=False, require_setup=True, de-
                                bug=False)
```

Saves the dict `data_dict` in the Photon-HDF5 format.

This function requires the data to be saved as `data_dict` argument. The data needs to have the hierarchical structure of a Photon-HDF5 file. For the purpose, we use a standard python dictionary: each key is a Photon-HDF5 field name and each value contains data (e.g. array, string, etc..) or another dictionary (in which case, it represents an HDF5 sub-group). Similarly, sub-dictionaries contain data or other dictionaries, as needed to represent the hierarchy of Photon-HDF5 files.

Features of this function:

- Checks that all field names are valid Photon-HDF5 field names.
- Checks that all field type match the Photon-HDF5 specs (scalar, array, or string).

- Populates automatically the identity group with filename, software, version and file creation date.
- Populates automatically the provenance group with info on the original data file (if it can be found on disk): creation and modification date, path.
- Computes field *acquisition_duration* when not provided (single-spot data only).

Minimal fields required to create a Photon-HDF5 file:

- /description* (string)
- /photon_data/timestamps* (array)
- /photon_data/timestamps_specs/timestamps_unit* (scalar float)
- /setup/num_pixels* (int): number of detectors
- /setup/num_spots* (int): number of excitation/detection spots
- /setup/num_spectral_ch* (int): number of detection spectral bands
- /setup/num_polarization_ch* (int): number of detected polarization states
- /setup/num_split_ch* (int): number of beam splitted channels
- /setup/modulated_excitation* (bool): True if excitation is alternated.
- /setup/lifetime* (bool): True if dataset contains TCSPC data.

See also [Writing Photon-HDF5 files](#).

As a side effect *data_dict* is modified by adding the key ‘_data_file’ containing a reference to the pytables file.

Parameters

- **data_dict** (*dict*) – the dictionary containing the photon data. The keys must strings matching valid Photon-HDF5 paths. The values must be scalars, arrays, strings or another dict.
- **h5_fname** (*string or None*) – file name for the output Photon-HDF5 file. If None, the file name is taken from *data_dict['_filename']* with extension changed to ‘.hdf5’.
- **user_descr** (*dict or None*) – dictionary of descriptions (strings) for user-defined fields. The keys must be strings representing the full HDF5 path of each field. The values must be binary (i.e. encoded) strings restricted to the ASCII set.
- **overwrite** (*bool*) – if True, a pre-existing HDF5 file with same name is overwritten. If False, save the new file by adding the suffix “new_copy” (and if a “_new_copy” file is already present overwrites it).
- **compression** (*dict*) – a dictionary containing the compression type and level. Passed to *pytables.tables.Filters()*.
- **close** (*bool*) – If True (default) the HDF5 file is closed before returning. If False the file is left open.
- **validate** (*bool*) – if True, after saving perform a validation step raising an error if the specs are not followed.
- **warnings** (*bool*) – if True, print warnings for important optional fields that are missing. If False, don’t print warnings.
- **skip_measurement_specs** (*bool*) – if True don’t print any warning for missing measurement_specs group.

- **require_setup** (*bool*) – if True, raises an error if some mandatory fields in /setup are missing. If False, allows missing setup fields (or missing setup altogether). Use False when saving only detectors' dark counts.
- **debug** (*bool*) – if True prints additional debug information.

For description and specs of the Photon-HDF5 format see: <http://photon-hdf5.readthedocs.org/>

```
phconvert.hdf5.assert_valid_photon_hdf5(datafile, warnings=True, verbose=False,
                                         strict_description=True, require_setup=True,
                                         skip_measurement_specs=False)
```

Asserts that `datafile` follows the Photon-HDF5 specs.

If the input datafile does not follow the specifications, it raises the `Invalid_PhotonHDF5` exception, with a message indicating the cause of the error.

This function checks that:

- all fields are valid Photon-HDF5 names
- all fields have valid descriptions
- all mandatory fields are present
- if /setup/lifetime is True (i.e. 1), assures that nanotimes and nanotimes_specs are present

Parameters

- **datafile** (*string* or *tables.File*) – input data file to be validated
- **warnings** (*bool*) – if True, print warnings for important optional fields that are missing. If False, don't print warnings.
- **verbose** (*bool*) – if True print details about the performed tests.
- **strict_description** (*bool*) – if True consider a non-conforming description (TITLE) a specs violation.
- **require_setup** (*bool*) – if True, raises an error if some mandatory fields in /setup are missing. If False, allows missing setup fields (or missing setup altogether).
- **skip_measurement_specs** (*bool*) – if True don't print any warning for missing measurement_specs group.

Utility functions

Utility functions to work with HDF5 files in pytables.

```
phconvert.hdf5.print_children(group)
```

Print all the sub-groups in *group* and leaf-nodes children of *group*.

Parameters *group* (*pytables group*) – the group to be printed.

```
phconvert.hdf5.print_attrs(node, which='user')
```

Print the HDF5 attributes for *node_name*.

Parameters

- **node** (*pytables node*) – node whose attributes will be printed. Can be either a group or a leaf-node.
- **which** (*string*) – Valid values are 'user' for user-defined attributes, 'sys' for pytables-specific attributes and 'all' to print both groups of attributes. Default 'user'.

`phconvert.hdf5.dict_from_group(group, read=True)`
Return a dict with the content of a PyTables *group*.

`phconvert.hdf5.dict_to_group(group, dictionary)`
Save *dictionary* into HDF5 format in *group*.

CHAPTER 2

Module loader

This module contains functions to load each supported data format. Each loader function loads data from a third-party formats into a python dictionary which has the structure of a Photon-HDF5 file. These dictionaries can be passed to `phconvert.hdf5.save_photon_hdf5()` to save the data in Photon-HDF5 format.

The loader module contains high-level functions which “fill” the dictionary with the appropriate arrays. The actual decoding of the input binary files is performed by low-level functions in other modules (`smreader.py`, `pqreader.py`, `bhreader.py`). When trying to decode a new file format, these modules can provide useful examples.

```
phconvert.loader.nsalex_bh(filename_spc, donor=4, acceptor=6, alex_period_donor=(10, 1500),
                           alex_period_acceptor=(2000, 3500), excitation_wavelengths=(5.32e-
07, 6.35e-07), detection_wavelengths=(5.8e-07, 6.8e-07), al-
low_missing_set=False, tcspc_num_bins=None, tcspc_unit=None)
```

Load a .spc and (optionally) .set files for ns-ALEX and return 2 dict.

The first dictionary can be passed to the `phconvert.hdf5.save_photon_hdf5()` function to save the data in Photon-HDF5 format.

Returns the first contains the main photon data (timestamps, detectors, nanotime, ...); the second contains the raw data from the .set file (it can be saved in a user group in Photon-HDF5).

Return type Two dictionaries

```
phconvert.loader.nsalex_ht3(filename, donor=0, acceptor=1, alex_period_donor=(150,
1500), alex_period_acceptor=(1540, 3050),
excitation_wavelengths=(5.23e-07, 6.28e-07),
detection_wavelengths=(5.8e-07, 6.8e-07))
```

Load a .ht3 file containing ns-ALEX data and return a dict.

This dictionary can be passed to the `phconvert.hdf5.save_photon_hdf5()` function to save the data in Photon-HDF5 format.

```
phconvert.loader.usalex_sm(filename, donor=0, acceptor=1, alex_period=4000, alex_offset=750,
alex_period_donor=(2850, 580), alex_period_acceptor=(930,
2580), excitation_wavelengths=(5.32e-07, 6.35e-07),
detection_wavelengths=(5.8e-07, 6.8e-07), software='LabVIEW
Data Acquisition usALEX')
```

Load a .sm us-ALEX file and returns a dictionary.

This dictionary can be passed to the `phconvert.hdf5.save_photon_hdf5()` function to save the data in Photon-HDF5 format.

This module contains functions to load and decode files from PicoQuant hardware.

The primary exported functions are:

- `load_ht3()` which returns decoded timestamps, detectors, nanotimes and metadata from an HT3 file.
- `load_pt3()` which returns decoded timestamps, detectors, nanotimes and metadata from a PT3 file.

Other lower level functions are:

- `ht3_reader()` which loads metadata and raw t3 records from HT3 files
- `pt3_reader()` which loads metadata and raw t3 records from PT3 files
- `process_t3records()` which decodes the t3 records returning timestamps (after overflow correction), detectors and TCSPC nanotimes.

Note that the functions performing overflow/rollover correction can take advantage of `numba`, if installed, to significantly speed-up the processing.

List of functions

High-level functions to load and decode several PicoQuant file formats:

`phconvert.pqreader.load_ht3(filename, ovfunc=None)`

Load data from a PicoQuant .ht3 file.

Parameters

- **filename** (*string*) – the path of the HT3 file to be loaded.
- **ovfunc** (*function or None*) – function to use for overflow/rollover correction of timestamps. If `None`, it defaults to the fastest available implementation for the current machine.

Returns A tuple of timestamps, detectors, nanotimes (integer arrays) and a dictionary with metadata containing at least the keys ‘`timestamps_unit`’ and ‘`nanotimes_unit`’.

`phconvert.pqreader.load_pt3(filename, ovfunc=None)`

Load data from a PicoQuant .pt3 file.

Parameters

- **filename** (*string*) – the path of the PT3 file to be loaded.
- **ovfunc** (*function or None*) – function to use for overflow/rollover correction of timestamps. If None, it defaults to the fastest available implementation for the current machine.

Returns A tuple of timestamps, detectors, nanotimes (integer arrays) and a dictionary with metadata containing at least the keys ‘timestamps_unit’ and ‘nanotimes_unit’.

Low-level functions

These functions are the building blocks for loading and decoding the different files:

`phconvert.pqreader.ht3_reader(filename)`

Load raw t3 records and metadata from an HT3 file.

`phconvert.pqreader.pt3_reader(filename)`

Load raw t3 records and metadata from a PT3 file.

`phconvert.pqreader.process_t3records(t3records, time_bit=10, dtime_bit=15, ch_bit=6, special_bit=True, ovfunc=None)`

Extract the different fields from the raw t3records array (.ht3).

Returns 3 arrays representing detectors, timestamps and nanotimes.

This module contains functions to load and decode files from Becker & Hickl hardware.

The high-level function in this module are:

- `load_spc()` which loads and decoded the photon data from SPC files.
- `load_set()` which returns a dictionary of metadata from SET files.

Becker & Hickl SPC Format

The structure of the SPC format is here described.

SPC-600/630

SPC-600/630 files have a record of 48-bit (6 bytes) in little endian (<) format. The first 6 bytes of the file are an header containing the *timestamps_unit* (in 0.1ns units) in the two central bytes (i.e. bytes 2 and 3). In the following drawing each char represents 2 bits:

```

bit: 64      48      0
    0000 0000 XXXX XXXX XXXX XXXX XXXX XXXX
        '-----' '---' '---' '-----'
field names:      a      c      b      d

    0000 0000 XXXX XXXX XXXX XXXX XXXX XXXX
        '-----' '---' '---' '-----'
numpy dtype:      a      c      b      field0

macrotime = [ b ] [      a      ] (24 bit)
detector   = [ c ]                (8 bit)
nanotime   = [ d ]                (12 bit)

overflow bit: 13, bit_mask = 2^(13-1) = 4096

```

SPC-134/144/154/830

SPC-134/144/154/830 files have a record of 32-bits (4 bytes) in little endian (<) format. The first 4 bytes of the file are an header containing the *timestamps_unit* (in 0.1ns units) in first two bytes. In the following drawing each char represents 2 bits:

```
bit:          32          0
             XXXX XXXX XXXX XXXX
             ' ' ' ' ' ' ' '
field names:  a      b      c      d

             XXXX XXXX XXXX XXXX
             ' ' ' ' ' ' ' '
numpy dtype:  field1    field0

macrotime = [ d ]      (12 bit)
detector  = [ c ]      (4 bit)
nanotime  = [ b ]      (12 bit)
aux       = [ a ]      (4 bit)

aux = [invalid, overflow, gap, mark]

If overflow == 1 and invalid == 1 --> number of overflows = [ b ][ c ][ d ]
```

List of functions

High-level functions to load and decode Becker & Hickl SPC/SET pair of files:

`phconvert.bhreader.load_spc(fname, spc_model='SPC-630')`

Load data from Becker & Hickl SPC files.

Parameters `spc_model` (*string*) – name of the board model. Valid values are 'SPC-630', 'SPC-134', 'SPC-144', 'SPC-154' and 'SPC-830'.

Returns 3 numpy arrays (timestamps, detector, nanotime) and a float (timestamps_unit).

`phconvert.bhreader.load_set(fname_set)`

Return a dict with data from the Becker & Hickl .SET file.

Low-level functions

These functions are the building blocks for decoding Becker & Hickl files:

`phconvert.bhreader.bh_set_identification(fname_set)`

Return a dict containing the IDENTIFICATION section of .SET files.

The both keys and values are native strings (binary strings on py2 and unicode strings on py3).

`phconvert.bhreader.bh_set_sys_params(fname_set)`

Return a dict containing the SYS_PARAMS section of .SET files.

The keys are native strings (traditional strings on py2 and unicode strings on py3) while values are numerical type or byte strings.

`phconvert.bhreader.bh_decode(s)`

Replace code strings from .SET files with human readable label strings.

`phconvert.bhreader.bh_print_sys_params` (*sys_params*)

Print a summary of the Becker & Hickl system parameters (.SET file).

CHAPTER 5

Indices and tables

- `genindex`
- `modindex`
- `search`

p

`phconvert.bhreader`, [11](#)
`phconvert.hdf5`, [3](#)
`phconvert.loader`, [7](#)
`phconvert.pqreader`, [9](#)

A

`assert_valid_photon_hdf5()` (in module `phconvert.hdf5`), [5](#)

B

`bh_decode()` (in module `phconvert.bhreader`), [12](#)

`bh_print_sys_params()` (in module `phconvert.bhreader`), [12](#)

`bh_set_identification()` (in module `phconvert.bhreader`), [12](#)

`bh_set_sys_params()` (in module `phconvert.bhreader`), [12](#)

D

`dict_from_group()` (in module `phconvert.hdf5`), [5](#)

`dict_to_group()` (in module `phconvert.hdf5`), [6](#)

H

`ht3_reader()` (in module `phconvert.pqreader`), [10](#)

L

`load_ht3()` (in module `phconvert.pqreader`), [9](#)

`load_pt3()` (in module `phconvert.pqreader`), [9](#)

`load_set()` (in module `phconvert.bhreader`), [12](#)

`load_spc()` (in module `phconvert.bhreader`), [12](#)

N

`nsalex_bh()` (in module `phconvert.loader`), [7](#)

`nsalex_ht3()` (in module `phconvert.loader`), [7](#)

P

`phconvert.bhreader` (module), [11](#)

`phconvert.hdf5` (module), [3](#)

`phconvert.loader` (module), [7](#)

`phconvert.pqreader` (module), [9](#)

`print_attrs()` (in module `phconvert.hdf5`), [5](#)

`print_children()` (in module `phconvert.hdf5`), [5](#)

`process_t3records()` (in module `phconvert.pqreader`), [10](#)

`pt3_reader()` (in module `phconvert.pqreader`), [10](#)

S

`save_photon_hdf5()` (in module `phconvert.hdf5`), [3](#)

U

`usalex_sm()` (in module `phconvert.loader`), [7](#)